


Video Package Assembly Instructions:

The Video Package is something that is becoming more and more common on the microscopes as they are purchased. So below you will find a quick step by step to attaching the video package.

The way that the Video and Digital Package is assembled is:

- 1) First remove the binocular head on the microscope.

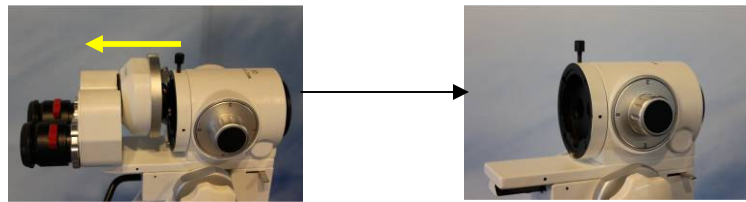

- 2) Second Attach the Beam Splitter:



- 3) Third insert the Video Adapter (20/80) and/or Digital Camera Adapter (50/50) into the Beamsplitter and tighten down the metal rings.

- 4) Attach Camera(s)

- 5) (Video) Attach TV to CCD Camera w/ S-video or composite cable

